

Instructions ActiveThree Biosemi AD-box

This room makes use of our newest Biosemi AD-box (ActiveThree). Please make sure to follow these instructions well as there are some small differences in the use.

1. On the desk outside the room, a battery charger is placed. To turn the ActiveThree AD-box on, you can grab ONE battery from the charging device and insert it at the side of the AD-box. Simply turn the knob, open the compartment, place the battery minus (-) first, then close the compartment. If you press the power button, it should turn on. (see pictures as a reference).
 - a. You can see how much the battery is charged on the charging device. If there are three solid green dots on top of a battery that is charging, it means it is full. If the green dots are blinking, it is still charging. The more dots, the more it is charged. A full charge should last anywhere from 10-20 hours.
 - b. Make sure to place the battery back into the charger and make sure it charges when you are finished!



2. The same electrode bundles and external electrodes can be used for the ActiveThree as the older ActiveTwo models. You can grab these from the usual place.
3. The CMS/DRL electrodes are **NOT** the same. You can find the new CMS/DRL electrodes hanging in the same place as before. There are two of them and each pair is labelled "Active Three Biosemi". They can be recognized by the red label and have a different input cable-ending (see pictures below).



PLEASE TURN OVER...

4. For the software, you will need to open an updated version of ActiView, specifically for the ActiveThree AD-boxes. On the PC, go to the file explorer and then: “*This PC*” → “*Local Disk (C:)*” → “*Biosemi*” → “*ActiView100-Beta3*”.
 - a. If you try to open the older version, it will not work!
5. Everything else is the same, good luck 😊