

# EyeLink Set-up

---

There are nine rooms with an EyeLink system available. The eye trackers have different models and specs. Please think of what you need for your experiment before you reserve a specific lab.


Example of one the Desktop Mount EyeLink cameras:



All nine EyeLink rooms have black stickers placed on the table. These stickers indicate where to place the equipment for the default set-up of: **eye-to-EyeLink distance of 55 cm**, and **eye-to-monitor distance of 70 cm**. If you prefer to use other distances, please move your equipment relative to these stickers without removing them. Please do not leave tape on the table to mark your personal preference. If you plan to use the same lab for a long period of time, please email [bblabs.fgb@vu.nl](mailto:bblabs.fgb@vu.nl) or the Lab Coordinator so we can place a second set of stickers on the table placed at your personal preference. When you are done using the lab, please place the equipment

back on the black stickers.

Visual representation of the default Eyelink set-up (table stickers):

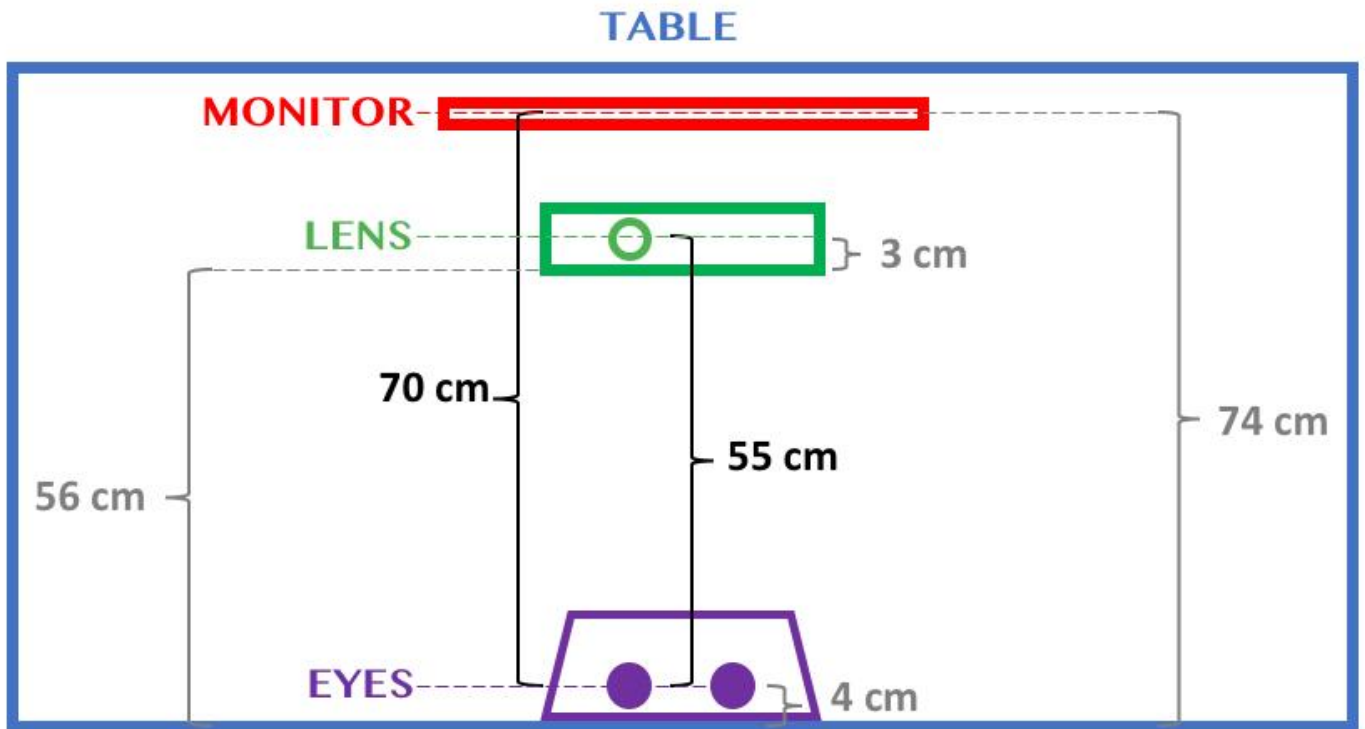

Revision #9

Created 2025-11-03 14:57:42 UTC by Admin

Updated 2026-06-08 13:43:17 UTC by Admin